

Touptek Microscope - HDMI-Camera Alpha-1080 HDMI + USB 2.0 + SD-Card Data transfer

ALPHA1080 series is a multiple outputs (HDMI+USB2.0+SD card) C-mount CMOS camera and it adopts ultra-high performance Sony CMOS sensor as the image-picking device. HDMI+USB2.0 are used as the data transfer interface to HDMI display or computer.

For HDMI camera mode, The XCamView will be loaded and the Camera Control Panel+ Measurement Toolbar and Synthesis Camera Control Toolbar are overlaid on the HDMI screen with the mouse move to the related region, in this case, the USB mouse can be used to set the camera, browse and compare the captured image, play the video and perform the measurement ital.

For USB Video camera mode, plug in the micro USB host cable to the camera's USB Video port and computer USB port, then the video stream can be transfer to computer with the advanced software ToupView/ToupLite. With ToupView/ToupLite, you can control the camera, process the video and image as ToupTek's other USB series camera.



The ALPHA1080 series' basic characteristics are as follows:

For HDMI output:

- All in 1(HDMI+USB+SD card) C-mount camera with Sony high sensitivity CMOS sensor
- Simultaneous HDMI & USB output
- Built-in mouse control
- Built-in image capture & video record to SD card
- Built-in camera control panel, including exposure(manual/auto)/gain, white balance(lockable), color adjustment, sharpness control
- Built-in video and image measurement
- Built-in toolbar including zoom, mirror, comparison, freeze, cross, browser functions
- Built-in image & video browsing, display & play
- Real time clock(RTC)

For USB Video mode:

- Ultra-Fine color engine with perfect color reproduction capability(USB)
- With advanced video & image processing application ToupView, which including professional image processing such as 2D measurement, HDR, image stitching, EDF(Extended Depth of Focus), image segmentation & count, image stacking, color composite and denoising(USB)
- Support standard UVC protocol for Windows/Linux/Mac(USB);
- CNC precision machining

The possible applications of ALPHA1080 series are as follows:

- Scientific research, education (teaching, demonstration and academic exchanges)
- Digital laboratory, medical research
- Industrial visual (PCB examination, IC quality control, mechanical part measurement)
- Medical treatment (pathological observation)
- Food (microbial colony observation and counting)
- Aerospace, military (high sophisticated weapons);




Order Code	Sensor & Size(mm)	Pixel(μm)	G Sensitivity Dark Signal	FPS/Resolution	Binning	Exposure
ALPHA1080A AL1080A	1080P/Sony IMX290 1/2.8"(5.57x4.13)	2.9x2.9	1300mv with 1/30s 0.15 mv with 1/30s	26@1920*1080P(HDMI) 26@1920*1080(USB)	1x1	0.036ms~8s
ALPHA1080B AL1080B	1080P/Sony IMX185 1/1.9"(7.20x4.50)	3.75x3.75	1120mv with 1/30s 0.15 mv with 1/30s	26@1920*1080P(HDMI) 26@1920*1080(USB)	1x1	0.34ms~4s

C: Color; M: Monochrome;

Interface & Button Functions		
	HDMI	HDMI Output
	USB Mouse	USB Mouse for the Control of the HDMI Mode
	USB Video	USB Video for the PC
	LED	Camera Status Indicate
	SD	SD Card Slot
	DC12V	12V/1A Power in
	ON/OFF	Power ON/OFF Switch
Other Specification for HDMI Output		
UI Operation	With USB Mouse for HDMI Mode	
Image Capture	High Speed in SD Card(8G)	
Video Record	1080P 26fps in SD Card(8G)(AVI Format in SD Card)	
Camera Control Panel	Including Exposure, Gain, White Balance, Color Adjustment and Sharpness Control	
Toolbar	Including Zoom, Mirror, Comparison, Freeze, Cross, Browser Function	
Other Specification for USB Output		
White Balance	Auto White Balance, Manual White Balance, ROI White Balance (👉)	
Color Technique	Ultra-Fine™ Color Engine (USB2.0)	
Capture/Control SDK	Standard UVC for Windows(USB2.0)	
Recording System	Still Picture or Movie (HDMI or USB)	
Software Environment (for USB2.0 Connection)		
Operating System	Microsoft® Windows® XP / Vista / 7 / 8 / 8.1 / 10(32 & 64 bit) OSx(Mac OS X) Linux	
PC Requirements	CPU: Equal to Intel Core2 2.8GHz or Higher	
	Memory:4GB or More	
	USB Port:USB2.0 High-speed Port	
	Display:19" or Larger CD-ROM	
Operating Environment		
Operating Temperature(in Centidegree)	-10~ 50	
Storage Temperature(in Centidegree)	-20~ 60	
Operating Humidity	30~80%RH	
Storage Humidity	10~60%RH	
Power Supply	DC 12V/1A Adapter	

Sizes of Alpha-1080

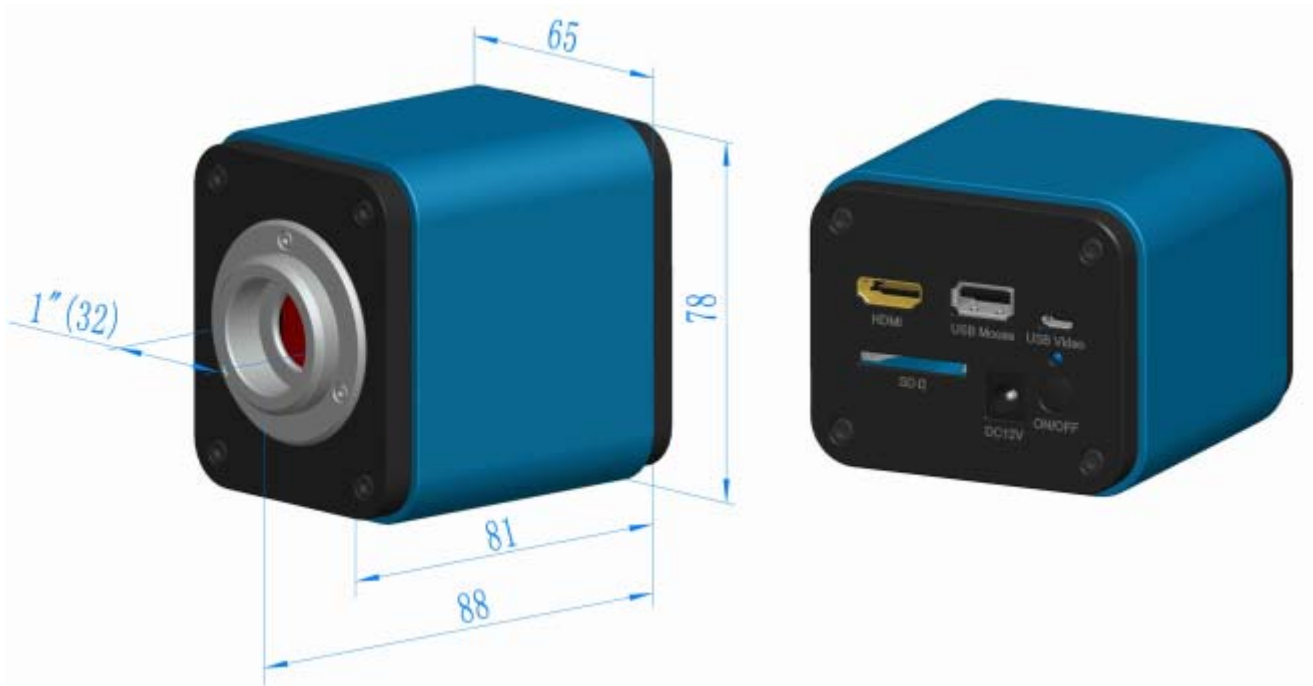

Product Images



ALPHA1080 Series and Its Back Panel




Different Views of ALPHA1080 Series



Mounted on microscope



With microscope and monitor

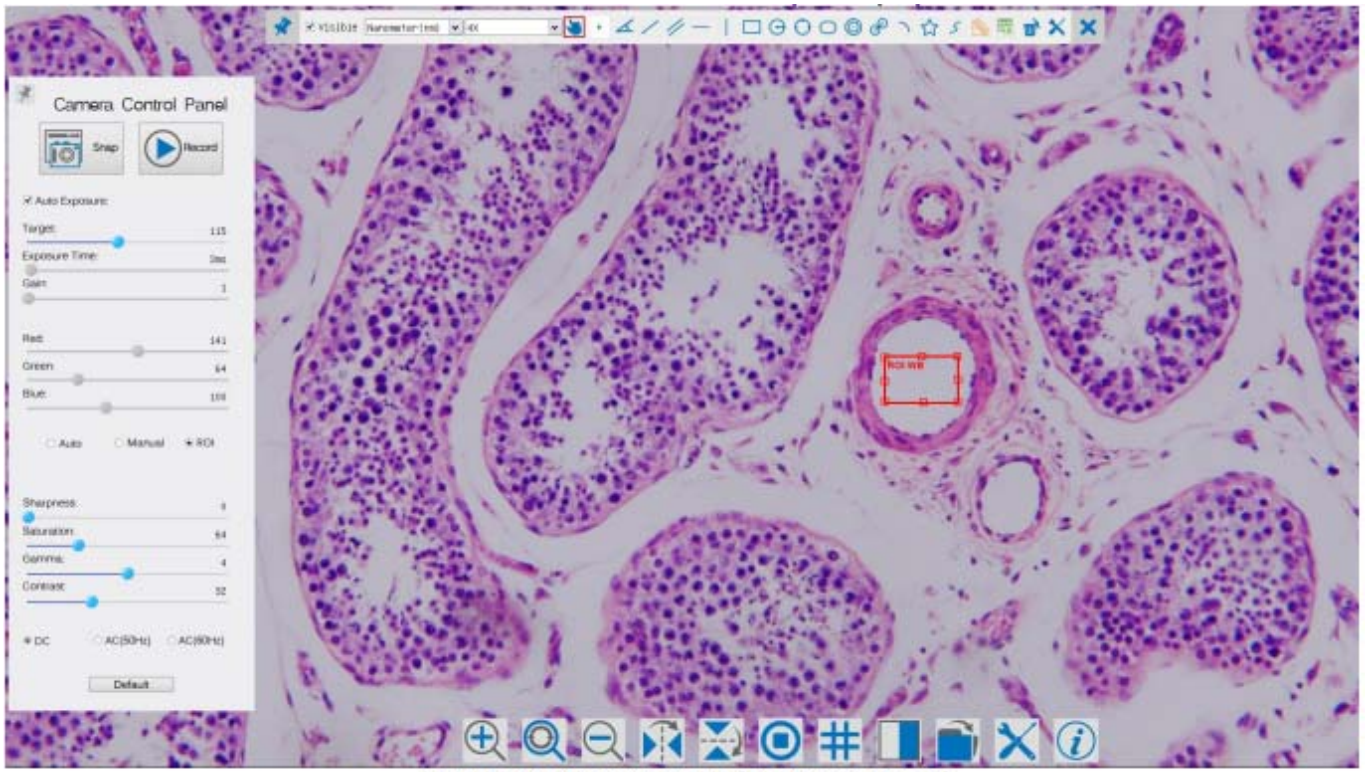

ALPHA1080 XCamView UI for Mouse Control

Packinmg information Alpha-1080



Adapters, glas scales, SD-card are not in scope of delivery

Standard Packing List			
A	Gift box : L:25.5cm W:17.0cm H:19.0cm (1pcs, 1.55kg/ box)		
B	ALPHA1080 Series camera		
C	Power Adapter: Input: AC 100~240V 50Hz/60Hz, Output: DC 12V 1A American standard: Model: GS12U12-P11 12W/12V/1A; UL/CUL/BSMI/CB/FCC EMI Standard:EN55022,EN61204-3, EN61000-3-2,-3, FCC Part 152 class B, BSMI CNS14338 EMS Standard: EN61000-4-2,3,4,5,6,8,11,EN61204-3,Class A Light Industry Standard European standard: Model:GS12E12-P11 12W/12V/1A; TUV(GS)/CB/CE/ROHS EMI Standard:EN55022,EN61204-3, EN61000-3-2,-3, FCC Part 152 class B, BSMI CNS14338 EMS Standard: EN61000-4-2,3,4,5,6,8,11,EN61204-3,Class A Light Industry Standard		
D	HDMI Cable		
E	USB Mouse		
F	Micro USB2.0 cable /2.0m		
G	CD (Driver & utilities software, Ø12cm)		
Optional Accessory			
H	Adjustable lens adapter	C-Mount to Dia.23.2mm eyepiece tube (Please choose 1 of them for your microscope)	108001/AMA037 108002/AMA050 108003/AMA075
		C-Mount to Dia.31.75mm eyepiece tube (Please choose 1 of them for your telescope)	108008/ATA037 108009/ATA050 108010/ATA075
I	Fixed lens adapter	C-Mount to Dia.23.2mm eyepiece tube (Please choose 1 of them for your microscope)	108005/FMA037 108006/FMA050 108007/FMA075
		C-Mount to Dia.31.75mm eyepiece tube (Please choose 1 of them for your telescope)	108011/FTA037 108012/FTA050 108013/FTA075
Note: For H and I optional items, please specify your camera type(C-mount, microscope camera or telescope camera), Touptek engineer will help you to determine the right microscope or telescope camera adapter for your application;			
J	108015(Dia.23.2mm to 30.0mm Ring)/Adapter rings for 30mm eyepiece tube		
K	108016(Dia.23.2mm to 30.5mm Ring)/ Adapter rings for 30.5mm eyepiece tube		
L	Calibration kit	106011/TS-M1(X=0.01mm/100Div.); 106012/TS-M2(X,Y=0.01mm/100Div.); 106013/TS-M7(X=0.01mm/100Div., 0.10mm/100Div.)	
M	SD Card(4G,8G,16G)		

The above statements are based on our present knowledge. Our statements should not be interpreted as a guarantee of characteristics. The use of our products by our customers is subject to different conditions, therefore none of our customers are relieved of the responsibility of testing our products by themselves. A liability for consequential damage will not be accepted in any case. For damage resulting from the use of this information we can only be held responsible if there is evidence of malice or negligence on our part. This data-sheet replaces any previous data sheets.

ASMETEC, METODRILL, METOCHECK, METOLIGHT, METOSTAT, METOCLEAN and METO are registered trade marks of ASMETEC GmbH.
 USB-CAM-ALPHA1080-DB-E.DOCX, VERSION MRZ-21